



YUCEBAS MACHINE ANALYTICAL EQUIPMENT INDUSTRY

## SEDIMENTATION DEVICE - Y15



YUCEBAS MACHINE ANALYTICAL EQUIPMENT INDUSTRY



### **SEDIMENTATION DEVICE - Y15**

SEDIMENTATION DEVICE determines the wheat bug damage in the wheat and flour samples. If there is a bug damage in the wheat, the sedimentation values will be low and it means lower protein quality. Sedimentation Device works for 5 minutes with 40 times shaking per minute with 30 degree angle on horizontal line according to the analysis principle. The users can do the tests with 6 different samples at the same time. It has a micro processor control and touchscreen LCD. The device can directly control from touch screen LCD. The device is suitable for Sedimentation Test and Modified Sedimentation Test. The users can do the Sedimentation Tests while waiting for the long resting period of the Modified Sedimentation Test. The internal temperature of the device, shaking number, test time and resting period of Modified Sedimentation Test as minutes while the users can do another Sedimentation Test, can shown on the touchscreen LCD.

#### **Technical Specifications :**

Dimensions : (H x D x W) 350 mm x 250 mm x 440 mm

Energy : 220 V 50 - 60 Hz

Net Weight : 14 Kg

### **YUCEBAS MACHINE ANALYTICAL EQUIPMENTS INDUSTRY**

**Factory :Fatih Mahallesi 451 Sokak No : 6**

**Helvacı / Aliğa / İZMİR / TURKEY**

**Tel : + 90 232 627 90 07 - Fax : + 90 232 627 90 08**

**[www.yucebasmakine.com.tr](http://www.yucebasmakine.com.tr)**

**[info@yucebasmakine.com.tr](mailto:info@yucebasmakine.com.tr)**

